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Description:

Quaternary ammonium compounds (QUATS) are often used in disinfectant solutions. To prepare the working solution, the disinfectant concentrate has to be diluted to the concentration for use. QUANTOFIX® QUAT facilitates checking and monitoring this concentration.

Pack content:

1 aluminum container with 100 test strips

Hazard warnings:

This test does not contain hazardous substances that must be labelled.

General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

Additional indications:

A greenish discoloration of the test strips when dry is normal and disappears after immersion in deionized water. It is recommended to assess the color of the test field in diffuse daylight.

Instructions for use:

1. Dip the test strip into the test solution for 5 seconds.
2. Shake off excess liquid.
3. Wait 15 seconds.
4. Compare with the color scale.

If QUATS are present, the test field turns green to bluish-green.

Note:

QUANTOFIX® QUAT can be used for the detection of quaternary ammonium compounds in general. The reference compound for the color scale is benzalkonium chloride. Other QUATS can be calculated as follows:

| Quaternary ammonium compound | Color scale value | | | | | | | | reaction time |
|--|--|---------------|---------------|---------------|----------------|----------------|----------------|-----------------|---------------|
| | 0 mg/L (ppm) | 10 mg/L (ppm) | 25 mg/L (ppm) | 50 mg/L (ppm) | 100 mg/L (ppm) | 250 mg/L (ppm) | 500 mg/L (ppm) | 1000 mg/L (ppm) | |
| | Corresponds to QUAT concentration (mg/L) | | | | | | | | |
| Benzalkonium chloride | 0 | 10 | 25 | 50 | 100 | 250 | 500 | 1000 | 15 sec |
| N-Cetyl-N,N,N-trimethylammonium bromide (CTAB) | 0 | 10 | 25 | 50 | 100 | 250 | 350 | 500 | 15 sec |
| N-cetylpyridinium bromide (= N-N-hexadecylpyridinium bromide) | 0 | 10 | 25 | 50 | 100 | 250 | 350 | 500 | 15 sec |
| Dodecyltrimethylammonium bromide (= Lauryltrimethylammonium bromide, LTAB) | 0 | 10 | 25 | 50 | 150 | 500 | 1500 | 2000 | 15 sec |
| Octadecyltrimethylammonium bromide | 0 | 10 | 25 | 50 | 100 | 250 | 500 | 1000 | 15 sec |
| Didecylmethylammonium bromide | 0 | 10 | 25 | 50 | 100 | 250 | 400 | 750 | 15 sec |
| Hyamine® 1622 ¹⁾ (= Benzethonium chloride) | 0 | 10 | 25 | 100 | 250 | 500 | 750 | 1500 | 60 sec |
| Cetyltrimethylammonium chloride (Benzyltrimethylhexadecylammonium chloride) | 0 | 10 | 25 | 50 | 100 | 250 | 500 | 1000 | 15 sec |
| Benzyltrimethylammonium bromide | 0 | 10 | 25 | 50 | 100 | 250 | 500 | 1000 | 60 sec |
| Dodecylguanidine monohydrochloride (N-2000 Antimicrobial) | 0 | 10 | 25 | 50 | 100 | 250 | 500 | 1000 | 15 sec |
| Preventol® DP1021 ²⁾ (= bronopol (INN), Dodecylguanidine monohydrochloride) | 0 | 5 | 10 | 25 | 50 | 100 | 250 | 500 | 15 sec |

¹⁾ Hyamine® is a registered trade mark of the Rohm and Haas Co

²⁾ Preventol® DP1021 is a registered trade mark of the Lanxess.

Conversion factor:

Concentration % \approx mg/L \div 10000

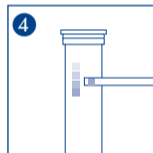
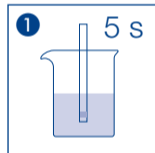
On request, we can produce our INDIQUAT test paper (REF 909000 – 909002) according to the customer's specifications.

Disposal:

Information regarding disposal can be found in the safety data sheet. You can download the SDS from www.mn-net.com/SDS.

Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 4 – 30 °C). If correctly stored, the test strips may be used until the use-by-date printed on the packaging.

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Schweiz: MACHEREY-NAGEL AG · Hirsackerstr. 7 · 4702 Oensingen · Schweiz